

LabVIEW™ Release Notes

LabVIEW 2010

These release notes contain installation instructions for LabVIEW and system requirements for the LabVIEW software.

If you are upgrading from a previous version of LabVIEW, read the *LabVIEW Upgrade Notes* included with the upgrade package before you install LabVIEW 2010. You must consider several issues before you convert VIs for use in LabVIEW 2010.

Read the *System Requirements* section of this document before you install LabVIEW and then follow the instructions in the *Installing LabVIEW 2010* section. After you install LabVIEW, read the *Where to Go from Here* section for information about getting started with and using LabVIEW.

Contents

System Requirements	2
Installing LabVIEW 2010.....	6
Windows	6
Mac OS X	7
Linux	8
Installing LabVIEW Add-Ons	9
Installing the Application Builder.....	9
Activating the LabVIEW License (Windows).....	10
Evaluating LabVIEW, Modules, or Toolkits.....	10
Single-Seat and Volume Licensing.....	11
Installing and Configuring Hardware	11
Windows	11
Mac OS X	11
Linux.....	12
Where to Go from Here	12
LabVIEW Quick Reference Card.....	12
ni.com	12

System Requirements

Table 1 describes the system requirements you need to run LabVIEW 2010.

Table 1. System Requirements for LabVIEW 2010

Platform	Media and System Requirements	Important Notes
All platforms	<p>LabVIEW requires a minimum of 256 MB of RAM, but National Instruments recommends 1 GB of RAM or higher.</p> <p>LabVIEW requires a screen resolution of at least 1,024 × 768 pixels.</p> <p>When you deploy a LabVIEW-built application, the LabVIEW Run-Time Engine requires a minimum of 64 MB of RAM and a screen resolution of 800 × 600 pixels for applications that require a screen, but National Instruments recommends 256 MB of RAM or higher and a screen resolution of 1,024 × 768 pixels or higher.</p>	<p>LabVIEW and the <i>LabVIEW Help</i> contain 16-bit color graphics. LabVIEW requires a minimum color palette setting of 16-bit color.</p> <p>You must have Adobe Reader installed to search PDF versions of all LabVIEW manuals.</p> <p>LabVIEW uses a directory for storing temporary files. National Instruments recommends that you have several megabytes of disk space available for this temporary directory. You can view or change the temporary directory by selecting Tools» Options and selecting Paths from the Category list. If LabVIEW aborts unexpectedly, it can leave files in the temporary directory. Remove old files from this directory to free disk space.</p> <p>(Windows) The default temporary directory is the system temporary directory, such as <code>C:\Documents and Settings\User\Local Settings\Temp</code>.</p> <p>(Mac OS X) Mac OS X 10.5.x uses <code>/private/var/folders/xxx/xxx/TemporaryItems</code>, where xxx is a secure, random name for each login.</p> <p>(Linux) The default temporary directory is <code>/tmp</code>.</p>

Table 1. System Requirements for LabVIEW 2010 (Continued)

Platform	Media and System Requirements	Important Notes
<p>Windows 7/ Vista/XP/ Windows Server 2003 R2 (32-bit)/ Windows Server 2008 R2 (64-bit)</p>	<p>LabVIEW requires a minimum of a Pentium III or greater or Celeron 866 MHz or equivalent processor, but National Instruments recommends at least a Pentium 4/M or equivalent processor. National Instruments recommends that you have at least 1.6 GB of disk space for the complete LabVIEW installation.</p> <p>When you deploy a LabVIEW-built application, the LabVIEW Run-Time Engine requires a minimum of a Pentium 200 MHz or equivalent processor, but National Instruments recommends a Pentium III or higher or Celeron 600 MHz or equivalent processor. The LabVIEW Run-Time Engine for controlling an application or front panel remotely requires at least 98 MB of disk space. The LabVIEW Run-Time Engine you include in a LabVIEW-built installer requires at least 217 MB of disk space. However, if you include any additional National Instruments installers in the LabVIEW-built installer you will require additional disk space.</p>	<p>LabVIEW does not support Windows 2000/NT/Me/98/95 or Windows XP x64.</p> <p>To use the <i>LabVIEW Help</i>, the Measurement & Automation Explorer (MAX) interactive help system, and the NI Example Finder, you must have Microsoft Internet Explorer 5.0 or later.</p> <p>To view and control a front panel remotely using Internet Explorer, National Instruments recommends that you have Internet Explorer 5.5 Service Pack 2 or later.</p>

Table 1. System Requirements for LabVIEW 2010 (Continued)

Platform	Media and System Requirements	Important Notes
Mac OS X	<p>LabVIEW requires Mac OS X 10.5.0 or later.</p> <p>This version of LabVIEW supports only Macintosh computers with Intel processors. LabVIEW does not support PowerPC processors. LabVIEW supports only Macintosh computers with case-insensitive HFS+ file systems.</p> <p>National Instruments recommends that you have at least 563 MB of disk space for the minimum LabVIEW installation or 1.2 GB of disk space for the complete LabVIEW installation.</p> <p>The LabVIEW Application Builder requires XCode 2.4.1 or later.</p>	<p>National Instruments recommends that you use Firefox 1.0.2 or later or Safari 1.3.2 or later to view the <i>LabVIEW Help</i>.</p>

Table 1. System Requirements for LabVIEW 2010 (Continued)

Platform	Media and System Requirements	Important Notes
Linux	<p>LabVIEW runs on Linux for Intel x86 processors with kernel version 2.2.x, 2.4.x, or 2.6.x.</p> <p>LabVIEW requires a minimum of a Pentium III or Celeron 866 MHz or equivalent processor, but National Instruments recommends a Pentium 4/M or equivalent processor.</p> <p>LabVIEW requires an X Window System server, such as XFree86 or X11R6.org.</p> <p>National Instruments recommends that you have at least 680 MB of disk space for the minimum LabVIEW installation, 890 MB of disk space for the complete LabVIEW installation, or 1.0 GB of disk space for the complete LabVIEW installation with drivers.</p> <p>LabVIEW is fully supported on Red Hat Enterprise Linux WS 4 or later and openSUSE 11.0 or later.</p> <p>LabVIEW can run without hardware driver support on any other distributions that provide GNU C Library (glibc, also known as libc.so.6) version 2.2.4 or later.</p>	<p>National Instruments recommends that you use Firefox 1.0.2 or later or Mozilla 1.2 or later to view the <i>LabVIEW Help</i>.</p> <p>LabVIEW requires GNU C Library version 2.2.4 or later. Most Linux vendors offer an updated <code>glibc rpm</code> for common Linux distributions on their Web sites.</p>

Installing LabVIEW 2010

Refer to the National Instruments Web site at ni.com/support if you encounter errors during installation.

Windows

The LabVIEW Platform DVDs include LabVIEW and the following modules and toolkits.

Data Management and Reporting

- Database Connectivity Toolkit
- DataFinder Toolkit
- Internet Toolkit
- Report Generation Toolkit for Microsoft Office

Embedded (Mobile and Robotics)

- Mobile Module
- Robotics Module

FPGA

- Compilation Tools for FPGA Devices other than Virtex-II
- Compilation Tools for Virtex-II FPGA Devices
- FPGA Compile Farm Toolkit
- FPGA Module

Image Acquisition and Image Processing

- Vision Development Module

Industrial Monitoring

- Datalogging and Supervisory Control Module
- Touch Panel Module

Interactive Datalogging and Measurements

- LabVIEW SignalExpress

Motion Control

- NI Motion Assistant
- NI SoftMotion Module

Real-Time

- Real-Time Execution Trace Toolkit
- Real-Time Module

Signal Processing, Analysis, and Math

- Adaptive Filter Toolkit
- Advanced Signal Processing Toolkit
- Digital Filter Design Toolkit
- MathScript RT Module
- Sound and Vibration

Simulation and Control

- Control Design and Simulation Module
- PID and Fuzzy Logic Toolkit
- Simulation Interface Toolkit
- System Identification Toolkit

Software Development and Deployment

- Desktop Execution Trace Toolkit
- Statechart Module
- Unit Test Framework Toolkit
- VI Analyzer Toolkit

If you purchased any of these add-ons, you can install them using the LabVIEW Platform DVDs. If you want to evaluate any of these add-ons before purchasing them, you can install these add-ons from the LabVIEW Platform DVDs and try them for free for 30 days.

Complete the following steps to install LabVIEW for Windows.

1. Log on as an administrator or as a user with administrator privileges.
2. Disable any automatic virus detection programs before you install. Some virus detection programs interfere with installation.
3. Insert the LabVIEW Platform DVD 1.



Note If you purchased your product with an NI Software Suite or NI Product Bundle, use the installation media that shipped with your purchase to install the product.

4. Follow the instructions on the screen to install and activate LabVIEW and any module or toolkit.



Note Some LabVIEW products have additional system requirements. As you select each product in the **Product List**, verify that the computer on which you are installing meets any additional requirements listed in the description field at the bottom of the window.

5. (Optional) Follow the instructions on the screen to install device drivers for National Instruments products. Refer to the *Installing and Configuring Hardware* section of this document for more information.
6. (Optional) Install any LabVIEW add-ons not found on the LabVIEW Platform DVDs. Refer to the *Installing LabVIEW Add-Ons* section of this document for more information.
7. After installation, enable any virus detection programs you disabled.
8. (Optional) Install and configure National Instruments hardware. Refer to the *Installing and Configuring Hardware* section of this document for more information.
9. Refer to the *Where to Go from Here* section of this document for more information about getting started with and using LabVIEW.



Note If you use Windows Vista or Windows XP, a **Security Alert** dialog box appears when you launch LabVIEW or the NI Example Finder for the first time. If you select the **Keep blocking this program** option, the LabVIEW VI Server, LabVIEW Web Server, and any server written in LabVIEW cannot accept incoming connections from a remote computer. Select the **Unblock this program, despite the security risk** option to configure your computer to launch LabVIEW without any changes in functionality. Refer to the National Instruments Web site at ni.com/info and enter the Info Code `expm69` for more information about correcting this problem.

To modify the current LabVIEW installation or to uninstall LabVIEW 2010, select **National Instruments Software** in the Add/Remove Programs applet in the Control Panel. When you modify the installation, a list of National Instruments software appears. Select a product in the list to add or remove individual components or to uninstall the product. Select multiple products to remove by pressing the <Shift> or <Ctrl> key while you select them. Click the **Uninstall** button to remove all the products you selected.

Mac OS X

Complete the following steps to install LabVIEW for Mac OS X.

1. Disable any automatic virus detection programs before you install. Some virus detection programs interfere with the installation program.
2. Insert the LabVIEW 2010 installation CD.
3. Run the appropriate installation program.

The following types of installations of LabVIEW 2010 are available:

- **Easy Install**—Installs all LabVIEW 2010 files, including LabVIEW, NI-488.2 drivers, and NI-VISA drivers. This is the default installation for LabVIEW.
 - **Custom Install**—If you select this option, you select the files to install. To select a custom installation, click the **Customize** button on the **Installation Type** page. You must select the **LabVIEW 2010** component to install the core set of LabVIEW files necessary for running LabVIEW 2010.
4. Follow the instructions that appear on the screen.
 5. After installation, enable any virus detection programs you disabled.
 6. (Optional) Install LabVIEW add-ons. Refer to the *Installing LabVIEW Add-Ons* section of this document for more information.

7. (Optional) Install and configure National Instruments hardware. Refer to the [Installing and Configuring Hardware](#) section of this document for more information.
8. Refer to the [Where to Go from Here](#) section of this document for more information about getting started with and using LabVIEW.

You cannot install LabVIEW into a non-default location, but you can move the `LabVIEW 2010` directory after you install. If you move the `LabVIEW 2010` directory to a non-default location, LabVIEW does not uninstall and remove the directory unless you move the `LabVIEW 2010` directory back to the default location.

To uninstall LabVIEW 2010, run the `Uninstall.sh` shell script file on the installation CD from a Terminal shell. The script prompts you to remove the files necessary to uninstall LabVIEW. If you want to keep any files you modified or added to the `LabVIEW 2010` directory, save those files in another location before uninstalling LabVIEW 2010.

Linux

Complete the following steps to install LabVIEW for Linux.

1. Log into the system as `root`.
2. Insert the LabVIEW 2010 installation CD. Use `mount /mnt/cdrom` to mount the CD. On some systems, the CD mounts automatically.
3. To change the current directory to the mounted CD, enter the following command:

```
cd /mnt/cdrom
```
4. To run the installation script, enter the following command:

```
sh ./INSTALL
```
5. (Optional) Install LabVIEW add-ons. Refer to the [Installing LabVIEW Add-Ons](#) section of this document for more information.
6. (Optional) Install and configure National Instruments hardware. Refer to the [Installing and Configuring Hardware](#) section of this document for more information.
7. Refer to the [Where to Go from Here](#) section of this document for more information about getting started with and using LabVIEW.



Note The format of the Academic Site License installation CD for Mac OS X and Linux is Mac Extended HFS+. If you are unable to mount the Academic Site License installation CD on a Linux system, use one of the following commands:

```
mount -t hfsplus /dev/scd0 /mnt/cdrom
mount -t hfsplus /dev/scd0 /media/cdrom
```

The `INSTALL` script installs LabVIEW in the `/usr/local` directory. The script also optionally runs the installation scripts for the NI-VISA and NI-488.2 drivers. The `INSTALL` script uses `rpm` to install on systems that support the `.rpm` format or to extract the `rpm` archives directly on other systems. You also can install the files manually by using `rpm`—or a GUI `rpm`-based install tool—on Red Hat or other `rpm`-based systems or on systems without `rpm` by using the utilities in the `bin` directory on the LabVIEW installation CD.

If you installed Netscape, Mozilla, or Firefox in the default location, the LabVIEW installation program automatically installs the browser plug-in so you can view and control front panels remotely using a Web browser. Otherwise, you can install the browser plug-in manually by installing the LabVIEW Run-Time Engine and then copying `/usr/local/lib/LabVIEW-2010/LV2010NPlugin.so` to the browser

plug-in directory such as `/usr/lib/netscape/plugins`, `/usr/lib/mozilla/plugins`, or `/usr/lib/firefox/plugins`.

Refer to the `README.html` or `README.txt` documents in `/mnt/cdrom` for custom installation instructions and other information.

Common LabVIEW Launch Errors on Linux

The following table lists common errors that might occur when you launch LabVIEW for Linux.

Error	Possible Cause and Solution
xlib: connection to :0.0 refused by server	Possible Cause —You are trying to run LabVIEW as a user who does not have permission to open a window on the display server. You typically receive these errors after you run the <code>su</code> command to temporarily become a different user, such as <code>root</code> (superuser). Solution —Exit the <code>su</code> command and launch LabVIEW as the login user or use the <code>xhost</code> or <code>xauth</code> commands to grant permission to open a window on the display server.
client is not authorized to connect to server	
internal error during connection authorization check	

Refer to the National Instruments Web site at ni.com/support for information about other errors that might occur when you launch LabVIEW for Linux.

Installing LabVIEW Add-Ons

Some versions of toolkits that are not found on the LabVIEW Platform DVDs do not work with LabVIEW 2010. Installing an incompatible toolkit might cause some features in the toolkit or LabVIEW to behave incorrectly. Installing modules and toolkits from the LabVIEW Platform DVDs ensures compatibility with LabVIEW 2010. Use the instructions in this section only to install add-ons not available on the LabVIEW Platform DVDs.

Complete the following steps to install LabVIEW add-ons.

1. Refer to the National Instruments Web site at ni.com/info and enter the Info Code `compat` to determine whether your LabVIEW add-on is compatible with the current version of LabVIEW.
2. Disable any automatic virus detection programs before you install. Some virus detection programs interfere with the installation program.
3. Insert the LabVIEW add-on installation CD and follow the instructions that appear on the screen.
4. After installation, enable any virus detection programs you disabled.

Refer to the National Instruments Web site at ni.com/labview/family for more information about National Instruments add-ons.

Installing the Application Builder

(Windows) LabVIEW installs the Application Builder when you install LabVIEW. If you activate the LabVIEW Professional Development System, you automatically activate the Application Builder. If you activate the LabVIEW Base Package or Full Development System, you must purchase and activate the Application Builder separately.

(Mac OS X and Linux) LabVIEW installs the Application Builder when you install the LabVIEW Professional Development System. If you install the LabVIEW Base Package or Full Development System, you must purchase the Application Builder separately.

Refer to the National Instruments Web site at ni.com/info and enter the Info Code `exd8yy` to access the Upgrade Advisor and purchase the Application Builder.

Activating the LabVIEW License (Windows)



Note (Mac OS X and Linux) LabVIEW does not require activation.

National Instruments uses activation to better support evaluation of our software, to enable additional software features, and to support license management in large organizations. To find out more about National Instruments software licensing, visit the National Instruments Web site at ni.com/license to find frequently asked questions, resources, and technical support.

You can activate licenses you have purchased for LabVIEW, modules, and toolkits during installation. If you do not activate a valid license for LabVIEW or any modules or toolkits that you install, the unlicensed product operates in evaluation mode. After you install the product, the evaluation begins the first time you launch the product and lasts 30 consecutive days. When the evaluation period expires, you must activate a valid license to continue using the product.

To activate the product license, use the serial number you received on the Certificate of Ownership included in the software kit. If your software kit does not include a Certificate of Ownership, you can find the serial number on the product packing slip or on the shipping label. If you have purchased the version of a product you want to activate, you also can use serial numbers from a previous version of the product to activate it. You can activate the LabVIEW license in any of the following ways:

- During installation, enter the serial number when prompted and complete the installation and activation process.
- After you launch LabVIEW in evaluation mode, click the **Activate Products** button in the **LabVIEW** dialog box.
- Select **Help»Activate LabVIEW Components** while LabVIEW is running in evaluation mode or select **Start»All Programs»National Instruments»NI License Manager** to launch NI License Manager. The license activation does not take effect until you restart LabVIEW.

If you do not activate LabVIEW, modules, or toolkits during installation, LabVIEW prompts you to activate when you launch LabVIEW. After you activate the licenses for all LabVIEW products you installed, you no longer see this prompt.

Evaluating LabVIEW, Modules, or Toolkits

You can access all the tools, VIs and functions, help, and examples available to the LabVIEW Professional Development System and most modules or toolkits in a free 30-day evaluation mode.

After the evaluation period for a module or toolkit expires, you are no longer able to perform the following actions:

- Run VIs from the module or toolkit.
- Run VIs that use VIs from the module or toolkit.
- Build VIs that use VIs from the module or toolkit into an executable or shared library.
- Access menus, tools, or palettes installed by the module or toolkit.

When you install a module or toolkit for evaluation, you install the help for the module or toolkit. The help for the module or toolkit remains installed after the 30-day evaluation period. Also, the NI Example Finder displays all module or toolkit examples after the 30-day evaluation period. You must uninstall the module or toolkit to remove the help and examples.

If you run LabVIEW in evaluation mode, you receive functionality of the Professional Development System. If you then purchase and activate a LabVIEW license, you receive full functionality of the development system you purchase. For example, if you create a VI using features that are available only in the Professional Development System and then you purchase and activate the Full Development System, the VI you created using the Professional Development System will be broken.

Single-Seat and Volume Licensing

LabVIEW supports both single-seat and volume licenses. Single-seat licensing is the use of LabVIEW by an individual user. Single-seat licenses may be used on one computer by many users or by one user on up to three computers. Volume licensing is available if you purchase five or more seats of the same product and requires a volume licensing agreement and additional volume licensing software. Each seat using LabVIEW must have a valid license from a license server. Refer to the National Instruments Web site at ni.com/license for more information about single-seat and volume licensing.

Installing and Configuring Hardware

After you install LabVIEW software, you can install the drivers you need to use National Instruments hardware. All National Instruments devices include the drivers you need to use the device. The drivers packaged with LabVIEW and those packaged with the devices might be different versions. LabVIEW works best with the latest available drivers that support the devices in your system. If you are using NI-DAQ 7.0 or later, refer to the *NI-DAQ Readme* for more information about which driver version to use with the data acquisition devices in your system.



Note National Instruments periodically updates device drivers. Refer to the National Instruments Web site at ni.com/info and enter the Info Code `exyckh` to download the latest drivers.

Install and configure National Instruments hardware after you install the drivers for those devices.

Windows

The NI Device Drivers media includes the drivers and other software you need to use all National Instruments hardware. Use MAX to confirm that LabVIEW recognizes the device, to configure accessories and device settings, and to run test panels to test device functionality, such as the ability to acquire and generate signals. Launch MAX by selecting **Start»All Programs»National Instruments»Measurement & Automation** or by double-clicking the **Measurement & Automation** icon on the desktop.

Refer to the installation instructions packaged with your device to install and configure your National Instruments device.

Mac OS X

The LabVIEW installation program installs the NI-VISA driver and the NI-488.2 driver for the GPIB interface by default.



Note (Mac OS X) LabVIEW does not currently support the full NI-DAQmx driver software. NI-DAQmx Base offers a subset of NI-DAQmx functionality on Mac OS X. Refer to the National Instruments Web Site at ni.com/info and enter the Info Code `exf8fb` to download NI-DAQmx Base.

Linux

The LabVIEW installation program prompts you to choose the NI-VISA and NI-488.2 drivers for the GPIB interface you are using. After you make a selection, LabVIEW installs those drivers automatically.



Note LabVIEW does not support the GPIB-1014 series (VME) devices or the original GPIB-SCSI box. LabVIEW does support the GPIB-SCSI-A box.

NI-DAQmx for Linux supports several National Instruments PXI, PCI, and SCXI data acquisition devices. Refer to the *NI-DAQmx Readme for Linux* for a complete list of supported devices. After you install the NI-DAQmx for Linux driver, refer to the *NI-DAQmx for Linux Configuration Guide* at `/usr/local/natinst/nidaqmx/docs/ConfigurationGuide.html` for information about testing and configuring a National Instruments DAQ device.

NI-DAQmx Base for Linux supports several National Instruments USB data acquisition devices. Refer to the *Readme for NI-DAQmx Base* for a complete list of supported devices. Documentation for the driver installs in the `/usr/local/natinst/nidaqmxbase/documentation` directory. Refer to the *LabVIEW Help* for more information about NI-DAQmx Base.

Where to Go from Here

Refer to the following resources for more information about LabVIEW, including new features and upgrade issues, programming concepts and step-by-step procedures, and known issues.

LabVIEW Quick Reference Card

The *LabVIEW Quick Reference Card* directs you to resources to reference during each phase of using LabVIEW—from installation through advanced development. This document references the *Getting Started with LabVIEW* manual, the *LabVIEW Upgrade Notes*, the *LabVIEW Help*, and other printed resources related to LabVIEW.

ni.com

Refer to ni.com/labviewzone for Web resources that include example VIs and discussion forums.

LabVIEW, National Instruments, NI, ni.com, the National Instruments corporate logo, and the Eagle logo are trademarks of National Instruments Corporation. Refer to the *Trademark Information* at ni.com/trademarks for other National Instruments trademarks. Other product and company names mentioned herein are trademarks or trade names of their respective companies. For patents covering National Instruments products/technology, refer to the appropriate location: **Help»Patents** in your software, the `patents.txt` file in your media, or the *National Instruments Patent Notice* at ni.com/patents