


# INSTALLATION GUIDE

## GPIB-USB-HS

To install the GPIB-USB-HS interface, connect the USB connector from the GPIB-USB-HS interface to an available USB port on your computer, as shown below. Before connecting the GPIB-USB-HS interface to GPIB devices, ensure that the computer and the GPIB devices are at the same ground potential<sup>1</sup>. The GPIB-USB-HS interface connects directly to most GPIB devices without requiring a GPIB cable. If your computer is already running, the operating system automatically detects the GPIB interface. Otherwise, the GPIB interface is detected when you start your computer.

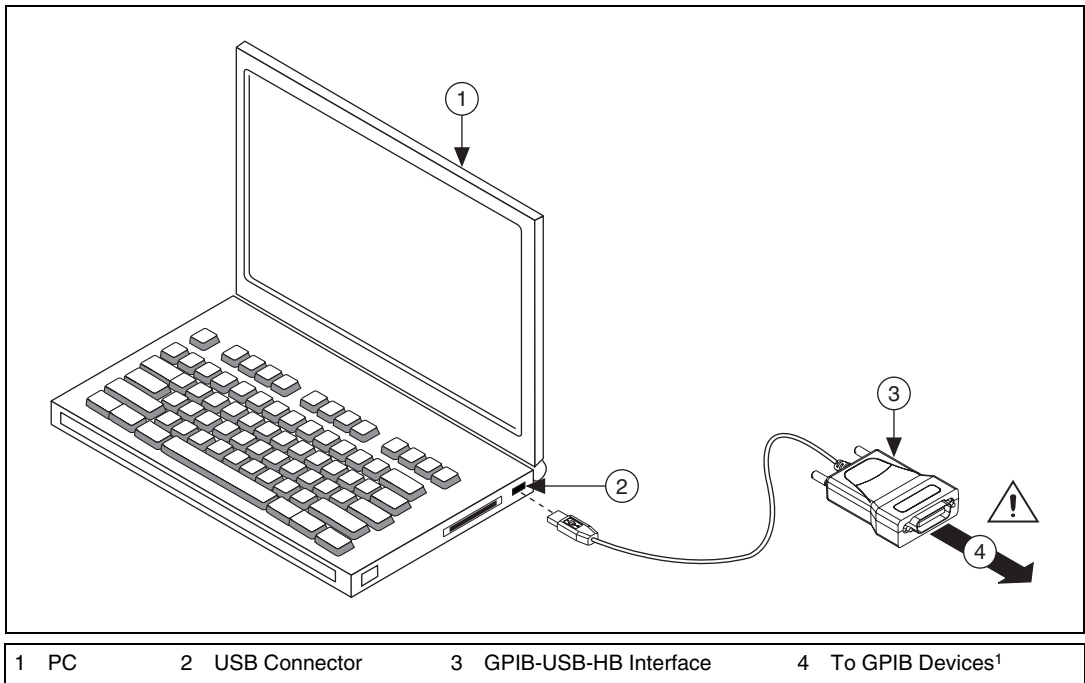

**Caution** <sup>1</sup>To connect to GPIB devices that do not share the same ground potential as the computer, use a GPIB bus expander/isolator or an isolated USB hub. Failure to isolate different ground potentials can damage the GPIB device, GPIB-USB interface, and computer.

National Instruments, NI, ni.com, and LabVIEW are trademarks of National Instruments Corporation. Refer to the *Terms of Use* section on [ni.com/legal](http://ni.com/legal) for more information about National Instruments trademarks. Other product and company names mentioned herein are trademarks or trade names of their respective companies. For patents covering National Instruments products/technology, refer to the appropriate location: **Help»Patents** in your software, the `patents.txt` file on your media, or the *National Instruments Patent Notice* at [ni.com/patents](http://ni.com/patents).